

3. -- -- Two dashes indicate the temperature, although still displayed, is above the range of the unit.

MOUNTING THE UNIT

1. To allow for different mounting positions, the white unit holder fits securely over the top or bottom of the unit.
2. Select the position and location where the unit will be mounted.
3. Make sure the surface of the location is clean and smooth.
4. Attach the unit to the selected location by removing the paper covering from the adhesive tape on the back of the unit holder and pressing the unit firmly in place.
5. The unit can be more permanently mounted by utilizing the two small screw holes on the unit holder. (Screws are not included.)

BENCH STAND

1. The bench stand is part of the BATTERY COMPARTMENT DOOR.
2. Locate the small rectangular opening at the bottom of the BATTERY COMPARTMENT DOOR.
3. Insert paper clip into opening and gently turn $\frac{1}{8}$ " in either direction. Stand will flip open.
4. To close, snap the stand closed.

ALL OPERATIONAL DIFFICULTIES

If this unit does not function properly for any reason, please replace the batteries with two new AAA alkaline batteries. (See "Battery Replacement" below.) Low battery power can occasionally cause any number of "apparent" operational difficulties. Replacing the batteries with fresh ones will solve most difficulties. Equivalent battery replacements are any AAA alkaline batteries.

BATTERY REPLACEMENT

1. To open battery compartment, slide BATTERY COMPARTMENT DOOR on the back of unit in direction of arrow.
2. Remove depleted batteries. Insert two fresh AAA alkaline batteries in appropriate positions to agree with the polarity symbols (+ and -) in the battery compartment.
3. Replace BATTERY COMPARTMENT DOOR.

TRACEABLE® HUMIDITY/ TEMPERATURE METER WITH DUAL MIN/MAX MEMORIES INSTRUCTIONS



WolfLabs

Pricing on any accessories shown can be found by keying the part number into the search box on our website.

The specifications listed in this brochure are subject to change by the manufacturer and therefore cannot be guaranteed to be correct. If there are aspects of the specification that must be guaranteed, please provide these to our sales team so that details can be confirmed.

www.wolflabs.co.uk

Tel : 01759 301142

Fax : 01759 301143

sales@wolflabs.co.uk

Please contact us if this literature doesn't answer all your questions.